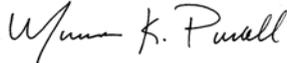


Standard Operating Procedure (SOP)

Title: Culture of *Renibacterium salmoninarum* from Tissues and Ovarian Fluid

Number: BACT-1

Version: 01 Created April 19, 2011

Approval:	Date:	Signature:
Bacteriology Supervisor:	April 30, 2011	

Effective Date:	April 30, 2011
Document Replaced:	
Reason for Revision:	
Area of Application:	

Purpose: Culture of *Renibacterium salmoninarum* in fish tissues or ovarian fluids

Sections:

- I. Important considerations before starting
- II. References
- III. Preparation of media and reagents
- IV. Culture procedure for kidney tissue
- V. Culture procedure for ovarian fluid

Disclaimer:

The use of trade, firm, or corporation names in this protocol is for the information and convenience of the reader. Such use does not constitute an official endorsement or approval by the U.S. Department of Interior or the U.S. Geological Survey of any product or service to the exclusion of others that may be suitable.

I. Important considerations before starting

Glassware: It is advisable to dedicate a separate set of glassware for culture to avoid antigen contamination of glassware used for other assays. This is most important in laboratories that also use immunological methods to detect *R. salmoninarum* antigen (e.g. ELISA). Extracellular proteins of the bacterium, especially the p57 antigen, can remain on container surfaces after conventional detergent washing and autoclaving.

pH meter: Check the pH of the medium before adding nurse medium to avoid contaminating probe with *R. salmoninarum* antigen.

II. References

- Evelyn, T.P.T. 1977. An improved growth medium for the kidney disease bacterium and some notes on using the medium. Bulletin of the Office of International Epizootics 87: 511-513.
- Evelyn, T.P.T., L. Prosperi-Porta, and J.E. Ketcheson. 1990. Two new techniques for obtaining consistent results when growing *Renibacterium salmoninarum* on KDM2 culture medium. Diseases of Aquatic Organisms 9: 209-212.
- Jansson, E., T. Hongslo, J. Höglund, and O. Ljungburg. 1996. Comparative evaluation of bacterial culture and two ELISA techniques for the detection of *Renibacterium salmoninarum* antigens in salmonid kidney tissues. Diseases of Aquatic Organisms 27:197-206.

III. Preparation of media and reagents

SKDM2: Recipe for 1 liter of selective kidney disease medium-2 (SKDM) agar plates (recipe makes approximately fifty 100 x 15 mm plates).

- Add measured amounts to a 2L flask
 - o Peptone 10 g
 - o Yeast extract 0.5 g
 - o Cysteine-HCL* 0.5 g
 - o Deionized water 900 mL
- Mix with stir bar and adjust medium to pH 6.5
- Add 15 mL of thawed nurse medium
- Add agar
 - o Agar 18g
- Autoclave medium on liquids cycle for 30-40 minutes at 15 psi and 121°C.
- Reduce temperature of the sterilized medium to 50°C
- Add 100 mL of fetal bovine serum (FBS)
- Add antibiotics (See Table 1: volume needed for 1 L)
- Dispense 20 mL agar per plate using a sterile disposable pipet
- Let plates cool for 15-20 minutes
- Seal plates into culture plate sleeve or plastic zipper bags and store in 15°C medium cabinet

Table 1: Antibiotics for SKDM2

Antibiotic	Stock solution ¹	Final w/v	Vol needed for 1 L
Cyclohexamide	1 g in 100 mL dH ₂ O ²	0.005%	5 mL
D-cycloserine	125 mg in 10 mL dH ₂ O	0.00125%	1 mL
Polymyxin-B sulfate	125 mg in 10 mL dH ₂ O	0.0025%	2 mL
Oxolinic acid	25 mg in 10 mL of 0.05N NaOH ³	0.00025	1 mL

¹Smaller volumes can be made. Antibiotics should be filter-sterilized (0.2 µm filter) before use.

²Cyclohexamide can be stored at 4°C but other antibiotics should be made fresh each time the medium is prepared.

³To make 0.05N NaOH: add 2.5 mL 1N NaOH in a 50 mL sterile volumetric flask and add sterile nH₂O to 50 mL mark.

*Note: The original protocol by Evelyn (1977) used 1 g/L cysteine. This concentration tends to result in crystal formation in the medium. Equivalent growth of *R. salmonarum* is attained *without* crystal formation if the cysteine is reduced to 0.5 g/L.

PBS-peptone: Recipe for 1 liter of 1X Voller phosphate buffered saline (PBS) pH 7.4 + 0.1% peptone.

- Add measured amounts into a 2 L bottle
 - o NaCl 8.0 g
 - o KH₂PO₄ 0.2 g
 - o Na₂HPO₄ anhydrous 1.09 g
 - o KCl 0.2 g
 - o Peptone 1.0 g
 - o Deionized H₂O to 1.0 L
- Mix with stir bar and check to make certain that pH is 7.4
- Autoclave on liquids cycle for 30-40 minutes at 15 psi and 121°C, then store in refrigerator at 4°C

Nurse Medium: Recipe for 1 liter of spent KDM-2 broth for use as a metabolite.

- Inoculate 1 liter of KDM 2 broth medium (recipe provided below) with 1 mL of stock *Renibacterium salmoninarum* isolate (ATCC 33209)
- Allow the culture to grow at 15° C for 10-14 days while being stirred
- Subject culture to centrifugation at 5000 x g for 20 minutes at 4°C
- Pool the supernatant and filter through a 0.2 µm bottle filter for sterilization.
 - o This may take several filters, depending on the amount of metabolite desired and how well the cells were separated out by centrifugation
- Dispense into 15 to 30-mL aliquots and freeze at -80° C

KDM2: Recipe for 1 liter of kidney disease medium-2 (KDM2) broth.

- Add measured amounts to a 2L flask
 - o Peptone 10 g
 - o Yeast extract: 0.5 g
 - o Cysteine-HCL* 0.5 g
 - o Deionized water 900 mL
- Mix with stir bar and adjust medium to pH 6.5
- Add thawed nurse medium
 - o Nurse medium 15 ml
- Autoclave medium on liquids cycle for 30-40 minutes at 15 psi and 121°C., then store in refrigerator at 4°C
- Reduce temperature of the sterilized medium to 50°C
- Add fetal bovine serum (FBS)
 - o FBS 100 mL
- Store in a 15°C medium cabinet or a 4°C refrigerator

*Note: The original protocol by Evelyn (1977) used 1 g/L cysteine. This concentration tends to result in crystal formation in the medium. Equivalent growth of *R. salmonarum* is attained *without* crystal formation if the cysteine is reduced to 0.5 g/L.

IV. Culture procedure for kidney tissue

Notes:

- *Keep tissue cold during entire procedure.*
- *Use freshly harvested kidney tissue.*

Method:

1. **Put the kidney sample into a pre-weighed sterile bag with a top closure.**
2. **Homogenize tissue with a heavy rolling pin or stomacher.** Homogenize small kidneys in a tube with a sterile pestle or other implement.
3. **Determine the tissue weight.**
Example: kidney tissue = 0.1 gram
4. **Add PBS-peptone at 10mL / gram tissue.** Mix tissue and PBS-peptone in bag by kneading the bag with fingers or by compression with a roller.
Example: 1ml for 0.1 gram tissue
5. **Transfer to a sterile centrifuge tube and subject to 2500 x g centrifugation for 20 minutes at 4°C.**
6. **Discard supernatant.**
7. **Resuspend pellet in PBS-peptone at 1:1 (w/v) ratio.** Vortex well.
Example: 0.1 gram of tissue would be resuspended in 100 µl (0.1 mL) PBS peptone
8. **Spread 10 µL of homogenate over entire surface SKDM2 agar plates with a sterile bent glass rod.** Alternatively, subject homogenates to serial 10-fold dilution in PBS-peptone for quantitative culture.
9. **Incubate at 15°C for up to 12 weeks.** Closure of plates with paraffin film or placement in a plastic zipper bag will retard drying.
10. **The presence of *R. salmoninarum* is typically confirmed by fluorescent antibody test (FAT) of colonies showing characteristic morphology.** See BACT-3 SOP for DFAT.
11. **Calculate CFU / g.** Divide the number of colonies by gram of tissue plated (take into account any further dilution for quantitative culture).
Example: If the starting tissue was 0.1 gram and 10 µl of homogenate was spread, than 0.01 g were used. So 10 colonies / 0.01 g = 1x10³ CFU/g

V. Procedure for Ovarian Fluid

Notes:

- *Keep ovarian fluid cold during entire procedure.*
- *Use freshly harvested ovarian fluid.*

Methods:

1. **Spread 100 µl of ovarian fluid over entire surface of SKDM2 agar plate with a sterile bent glass rod.** Alternatively, subject homogenates to serial 10-fold dilution in PBS-peptone for quantitative culture.
2. **Incubate at 15°C for up to 12 weeks.** Closure of plates with paraffin film or placement in a plastic zipper bag will retard drying.
3. **The presence of *R. salmoninarum* is typically confirmed by fluorescent antibody test (FAT) of colonies showing characteristic morphology.** See BACT-3 SOP for DFAT.
4. **Calculate CFU / mL.** Divide the number of colonies by mL of tissues plated (take into account any further dilution for quantitative culture).

Example:

If the starting volume was 0.1 mL and 100 colonies were observed:

$$100 \text{ colonies} / 0.1 \text{ mL} = 1 \times 10^3 \text{ CFU/mL}$$

If the ovarian fluid was diluted 1:10 before plating, then:

$$(100 \text{ colonies} / 0.1 \text{ mL}) * 10 \text{ fold dilution} = 1 \times 10^4 \text{ CFU/mL}$$